



抗体藥物複合体 (部位特異的ADC)
Antibody drug conjugate (site-specific ADC)

U241217A

Before drug conjugation

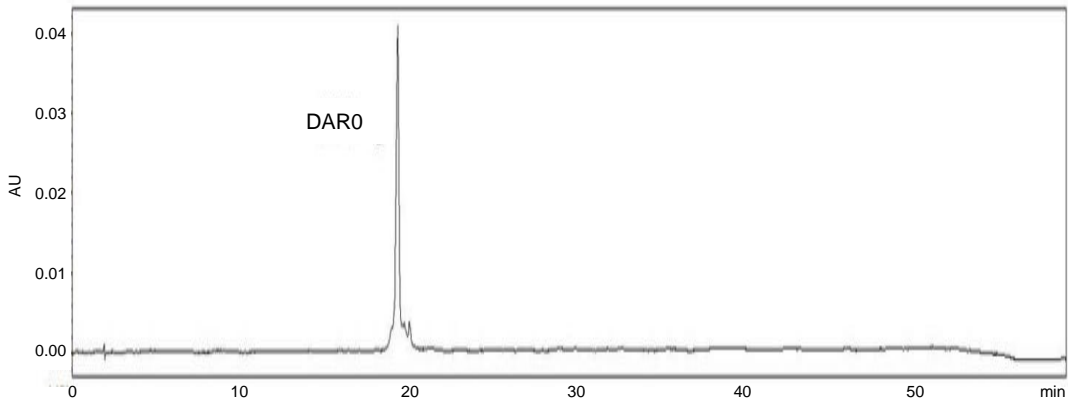

After drug conjugation

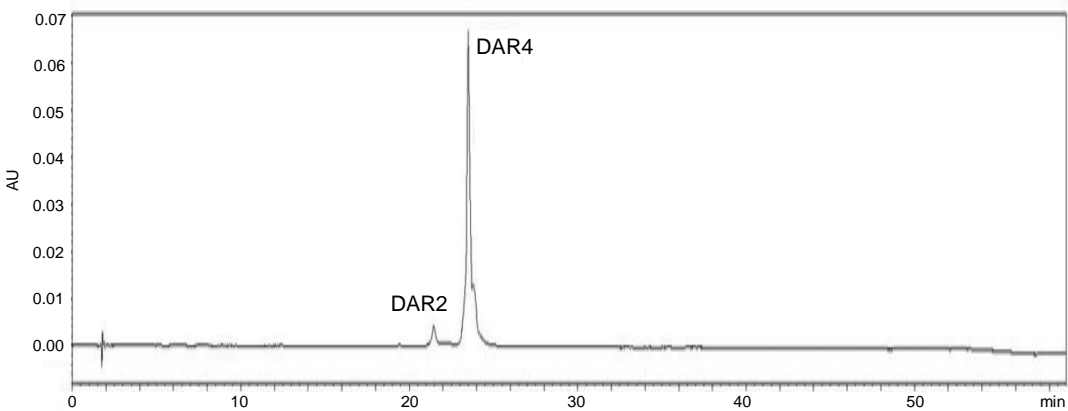

Courtesy of KriSan Biotech Co., Ltd.

Column : YMC-Triart Bio C4 (3 μ m, 30 nm)
150 X 4.6 mm I.D.

Eluent : A) water/TFA (100/0.1)
B) acetonitrile/TFA (100/0.1)
30%B (0–2.2 min), 30–46%B (2.2–44 min), 46–90%B (44–53 min), 90%B (53–59 min)

Flow rate : 1.0 mL/min

Temperature : 60°C

Detection : UV at 280 nm

Injection : 0.7 μ L (20 mg/mL)

Sample : Site-specific, 3rd generation ADC (in-house)