


ブテナフィン塩酸塩液・スプレー・クリーム


Butenafine hydrochloride solution and spray and cream

R100830A

Commercially available reagent
(0.5, 0.17 mg/mL, correspond to standard solution
for butenafine hydrochloride cream)

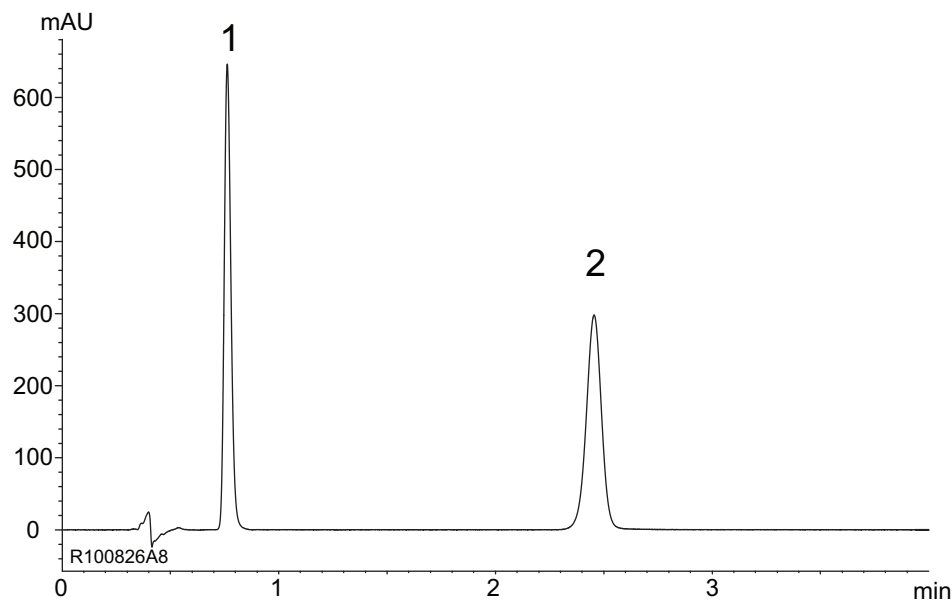


1.

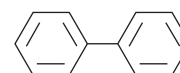


Diphenyl (I.S.)

2.

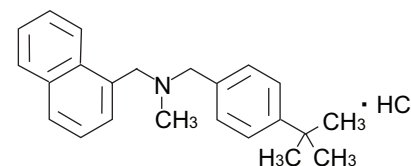

Butenafine hydrochloride

Sample solution for butenafine hydrochloride cream

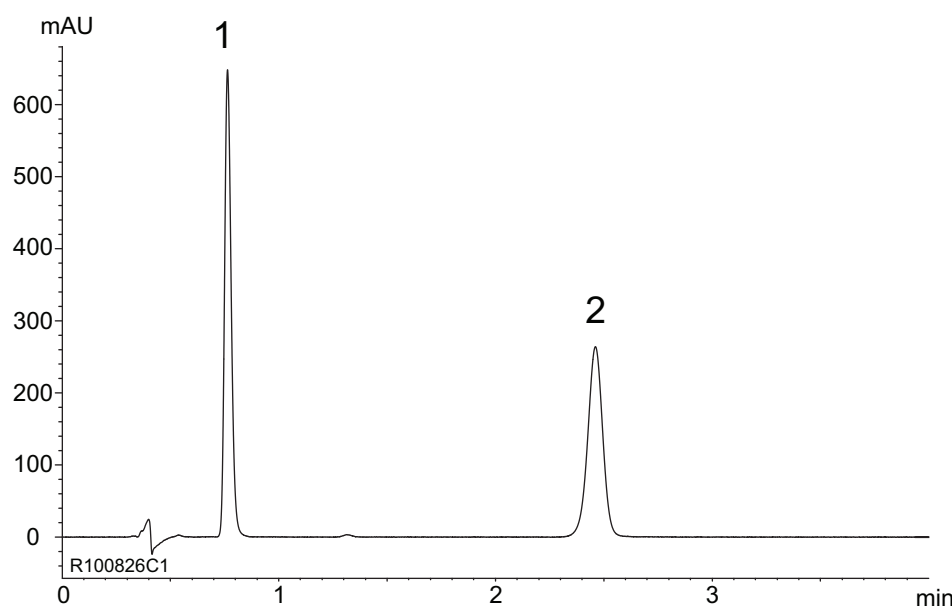

Column : YMC-Pack Pro C8 (3 μ m, 12 nm)
50 X 3.0 mm I.D.

Eluent : 1 mM CH₃COONH₄/acetonitrile (1/4)

Flow rate : 0.7 mL/min (adjust the flow rate so that the retention time of butenafine is about 2.5 min)

Temperature : 40°C

Detection : UV at 282 nm

Injection : 5 μ L

(日本薬局方収載原案記載条件)